



## Advanced flow cytometry

# FASER Kit – APC

Order No. 130-091-762

### Index

1. Description
  - 1.1 Background and product applications
  - 1.2 Reagent requirements
2. General protocol for immunofluorescent staining
3. Example of immunofluorescence amplification using the FASER Kit – APC

### 1. Description

<b>Components</b>	2×2 mL FcR Blocking Reagent, 1 mL APC-Activator (Reagent 1), 1 mL APC-Enhancer (Reagent 2).
<b>Size</b>	For 10 <sup>9</sup> total cells, up to 100 stainings.
<b>Staining concentration</b>	FcR Blocking Reagent is used at a dilution of 1:5. The APC-Activator (Reagent 1) and the APC-Enhancer (Reagent 2) are used at a dilution of 1:11.
<b>Product format</b>	Reagents are supplied as a solution containing 0.1% gelatine and 0.05% sodium azide.
<b>Storage</b>	Store protected from light at 4–8 °C. Do not freeze. The expiration date is indicated on the vial label.

#### 1.1 Background and product applications

The FASER (Fluorescence Amplification by Sequential Employment of Reagents) Kit – APC (allophycocyanin) is developed to increase the fluorescence intensity of cells labeled with an APC-conjugated antibody. The intensity of the APC fluorescence is amplified by sequential addition of the APC-Activator (Reagent 1) and the APC-Enhancer (Reagent 2). For further enhancement of the fluorescence intensity, sequential addition of the APC-Activator (Reagent 1) and the APC-Enhancer (Reagent 2) can be repeated several times.

The FASER Kit – APC does not affect direct staining with other fluorochrome-conjugated antibodies, but should not be combined with indirect immunofluorescent staining using biotin-conjugated antibodies. The Kit is suitable for APC-labeled, fresh or formaldehyde-fixed cells of any type or species in suspension. Analysis is performed by flow cytometry.


#### Examples of product application

- Amplification of APC fluorescence which is weak, for example
  - due to APC-conjugated antibody staining of antigens which are expressed at low levels,
  - due to immunomagnetic and APC-conjugated antibody labeling of the same antigen epitope,
  - due to APC-conjugated antibody staining with a low affinity antibody.
- Amplification of the fluorescence signal of an APC-labeled cell fraction for a clearer flow cytometric discrimination from the non-labeled cell fraction.
- Amplification of magnetic labeling with Anti-APC MicroBeads (# 130-090-855) by increasing the number of Anti-APC MicroBead binding sites per cell.

#### 1.2 Reagent requirements

- Buffer: PBS (phosphate buffered saline) pH 7.2, supplemented with 0.5% BSA (bovine serum albumin) and 2 mM EDTA. Keep buffer cold (4–8 °C).

▲ **Note:** EDTA can be replaced by other supplements such as anticoagulant citrate dextrose formula-A (ACD-A) or citrate phosphate dextrose (CPD). BSA can be replaced by other proteins such as serum albumin, serum or fetal calf serum. Buffers or media containing Ca<sup>2+</sup> or Mg<sup>2+</sup> are not recommended for use.

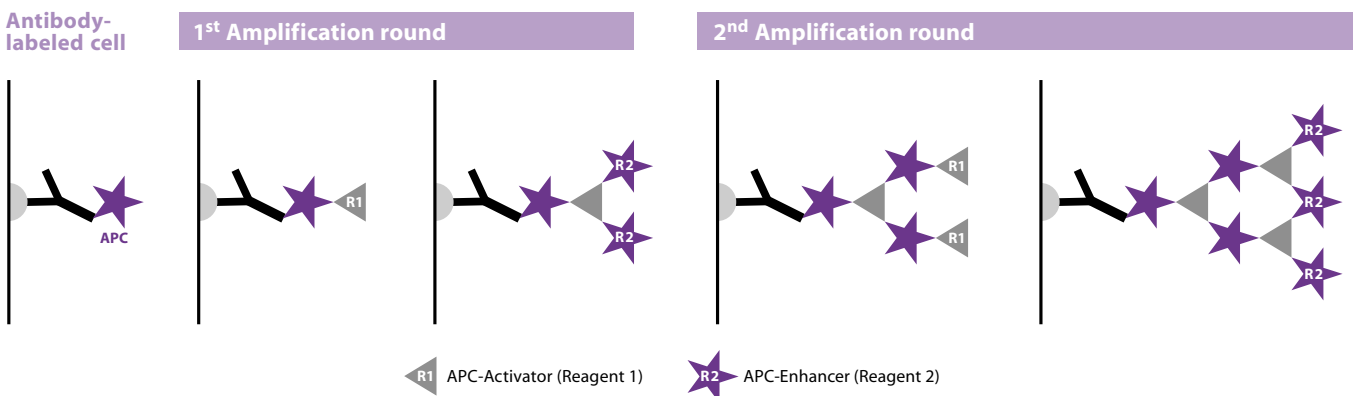

140-001-181-01

**Miltenyi Biotec**

www.miltenyibiotec.com



Miltenyi Biotec GmbH  
Friedrich-Ebert-Str. 68  
51429 Bergisch Gladbach, Germany  
Phone +49-2204-8306-0 Fax +49-2204-85197

Miltenyi Biotec Inc.  
12740 Earhart Avenue, Auburn CA 95602, USA  
Phone 800 FOR MACS, 530 888-8871  
Fax 530 888-8925

## 2. General protocol for immunofluorescent staining

▲ If formaldehyde fixation of cells is required, cells should be stained and treated with the FASER Kit – APC prior to fixation.


▲ If the FASER Kit – APC is used for amplification of magnetic labeling with Anti-APC MicroBeads, amplify APC staining as described below. For subsequent magnetic labeling, refer to the Anti-APC MicroBead data sheet.

▲ Volumes for fluorescent labeling given below are for up to  $10^7$  total cells. When working with fewer than  $10^7$  cells, use the same volumes as indicated. When working with higher cell numbers, scale up all reagent volumes and total volumes, accordingly (e.g. for  $2 \times 10^7$  total cells, use twice the volume of all indicated reagent volumes and total volumes).

1. Label cells with desired APC-conjugated antibody and afterwards wash cells according to the manufacturer's recommendations.
2. Resuspend up to  $10^7$  cells in 80  $\mu$ L of buffer.
3. Add 20  $\mu$ L of **FcR Blocking Reagent**.
4. Add 10  $\mu$ L of **APC-Activator (Reagent 1)**.
5. Mix well and incubate for 10 minutes in the dark at 4–8 °C.
  - ▲ **Note:** Working on ice requires increased incubation time. Higher temperatures and/or longer incubation times lead to non-specific cell labeling.
6. Wash cells by adding 1–2 mL of buffer per  $10^7$  cells and centrifuge at 300 $\times$ g for 10 minutes. Pipette off supernatant completely.
7. Resuspend cell pellet in 80  $\mu$ L of buffer.
8. Add 20  $\mu$ L of **FcR Blocking Reagent**.
9. Add 10  $\mu$ L of **APC-Enhancer (Reagent 2)**.
10. Mix well and incubate for 10 minutes in the dark at 4–8 °C.
  - ▲ **Note:** Working on ice requires increased incubation time. Higher temperatures and/or longer incubation times lead to non-specific cell labeling.
11. Wash cells by adding 1–2 mL of buffer per  $10^7$  cells and centrifuge at 300 $\times$ g for 10 minutes. Pipette off supernatant completely.
12. (Optional) Repeat steps 2–11, if further amplification of the APC fluorescence is required.
13. Finally resuspend cell pellet in a suitable amount of buffer for analysis by flow cytometry.

## 3. Example of immunofluorescence amplification using the FASER Kit – APC

Human peripheral blood mononuclear cells (PBMC) were stained with CD19-APC, human (# 130-091-248), and the APC fluorescence was increased by two rounds of amplification using the FASER Kit – APC. Cells were analyzed by flow cytometry. Cell debris and dead cells were excluded from the analysis based on scatter signals and propidium iodide (PI) fluorescence.

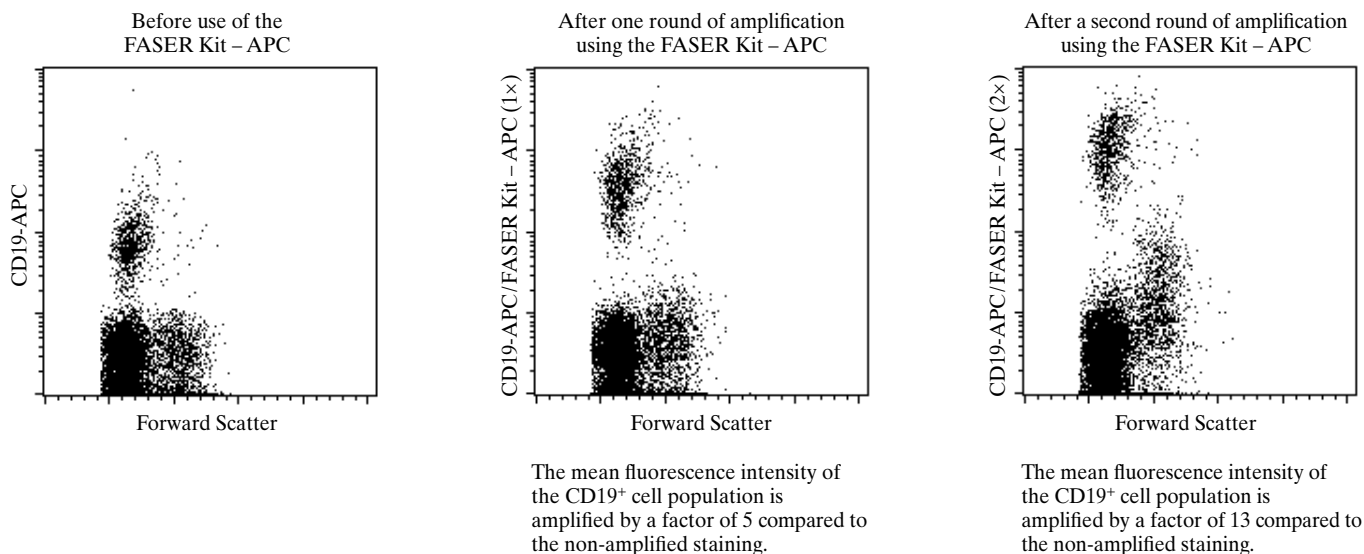

### Warning

Reagents contain sodium azide. Sodium azide yields hydrazoic acid under acid conditions, which is extremely toxic. Azide compounds should be diluted with running water before discarded. These precautions are recommended to avoid deposits in plumbing where explosive conditions may develop.

### Warranty

The products sold hereunder are warranted only to be free from defects in workmanship and material at the time of delivery to the customer. MILTENYI BIOTECH GmbH makes no warranty or representation, either expressed or

implied, with respect to the suitability of a product for a particular purpose. There are no warranties, expressed or implied, which extend beyond the Technical Specifications of the products. MILTENYI BIOTECH GmbH's liability is limited to either replacement of the products or refund of the purchase price.

MILTENYI BIOTECH GmbH is not liable for any property damage, personal injury or economic loss caused by the product.

MACS<sup>®</sup> is a registered trademark of Miltenyi Biotec GmbH.