

Contents

1. Description

1.1 Principle of the Epidermis Dissociation Kit ACF

1.2 Background information

1.3 Applications

1.4 Reagent and instrument requirements

2. Protocol

2.1 Reagent preparation

2.2 Epidermis dissociation protocol

2.2.1 Separation of epidermis from dermis (day 1)

2.2.2 Dissociation of epidermis (day 2)

2.2.3 Automated dissociation using the gentleMACS Dissociator

1. Description

| | |
|-------------------|--|
| Components | 4 vials, containing 13 mL of Buffer S (20× Stock Solution) 1 vial of Enzyme G, ACF (lyophilized powder) 2.5 mL of Enzyme P, ACF 1 vial of Enzyme A, ACF (lyophilized powder) |
| Size | For 25 digestions. The specified number of digestions is valid when digesting skin with an average weight of 1.5 g following the protocol in chapter 2.2. |
| Storage | Upon arrival immediately store Enzyme P, ACF in aliquots at -20 °C. Store all other components at 2–8 °C upon arrival. Reconstitute Enzymes G, ACF and Enzyme A, ACF before the date indicated on the box label. For information about reconstitution and storage after reconstitution of the lyophilized components refer to chapter 2.1. |

1.1 Principle of the Epidermis Dissociation Kit ACF

Epidermal tissue can be dissociated to single-cell suspensions by combining mechanical dissociation with enzymatic degradation of the extracellular adhesion proteins which maintain the structural integrity of tissues.

In a first step the epidermal cell layer is removed from the dermal cell layer after enzymatic treatment over night at 4 °C. The epidermal tissue is then further digested enzymatically and dissociated into a single-cell suspension by either using the gentleMACS™ Dissociators or by manual handling. Cells should be processed immediately for downstream applications, such as cell separation, cellular, or molecular analysis.

1.2 Background information

The Epidermis Dissociation Kit ACF, mouse enables the gentle and efficient generation of single-cell suspensions from mouse epidermal tissue. It has been particularly developed for the isolation of epidermal Langerhans cells.

This product is animal component-free (ACF) as it does not contain any primary raw materials derived directly from animals during the manufacturing process.

1.3 Applications

- Dissociation of mouse skin tissue for the isolation of Langerhans cells using the Epidermal Langerhans Cell MicroBead Kit, mouse (# 130-095-408)
- Phenotyping or enumeration of epidermal cell populations by flow cytometry

1.4 Reagent and instrument requirements

- Phosphate-buffered saline (PBS), pH 7.2
- PB buffer: Prepare a solution containing PBS, pH 7.2 and 0.5% bovine serum albumin (BSA) by diluting MACS® BSA Stock Solution (# 130-091-376) 1:20 with PBS. Keep buffer cold (2–8 °C).
- MACS SmartStrainers (70 µm) (50 filters, #130-098-462; 4×25 filters, #130-110-916)
- (Optional) gentleMACS Dissociator (# 130-093-235) or gentleMACS Octo Dissociator with Heaters (# 130-096-427)
- gentleMACS C Tubes (25 tubes, # 130-093-237; 4×25 tubes, # 130-096-334)

2. Protocol

▲ For details on the use of the gentleMACS Dissociators, refer to the gentleMACS Dissociator user manuals.

▲ For cell culture experiments subsequent to tissue dissociation, all steps should be performed under sterile conditions.

▲ The protocol is developed for the digestion of skin derived from 6–10 weeks old female C57 BL/6 or Balb/c mice.

▲ **Note:** It is not recommended to use skin derived from younger or older mice.

▲ An amount of up to 1.5 g skin tissue can be used per digestion, which corresponds approximately to the weight of back skin derived from two mice. When working with less than 1.5 g, use the same volumes as indicated below.

2.1 Reagent preparation

▲ Prepare 1× Buffer S by adding, for example, 1 mL of 20× Buffer S aseptically to 19 mL of sterile, distilled water. Store at 2–8 °C.

▲ **Note:** Handle under sterile conditions.

▲ Prepare Enzyme G, ACF by reconstitution of the lyophilized powder in the vial with 3 mL of sterile, distilled water. Do not vortex. Produce aliquots of appropriate volume. Store aliquots at –20 °C. Avoid repeated freeze-thaw-cycles.


▲ Prepare aliquots of appropriate volume of Enzyme P, ACF. Store aliquots at –20 °C. Avoid repeated freeze-thaw-cycles.

▲ Prepare Enzyme A, ACF by reconstitution of the lyophilized powder in the vial with 0.25 mL of sterile, distilled water. Do not vortex. Produce aliquots of appropriate volume. Store aliquots at –20 °C. Avoid repeated freeze-thaw-cycles.

2.2 Epidermis dissociation protocol


2.2.1 Separation of epidermis from dermis (day 1)

1. Remove hair completely from the back of the mice by plucking against the growth direction of the hair.
2. Cut skin at the base of the tail with a transverse cut using scissors and incise then upwards on both sides. Peel the tissue off with tweezers and make a transverse cut at the neck.
3. Wash tissue by placing the skin in a 100 mm petri dish containing 20 mL of PBS buffer.
4. Determine weight of the tissue to make sure that the limit of 1.5 g per digestion is not exceeded.
5. Transfer tissue to a new 100 mm petri dish with the epidermal side facing downwards.
6. Carefully scrape off the subcutaneous fat using a scalpel.

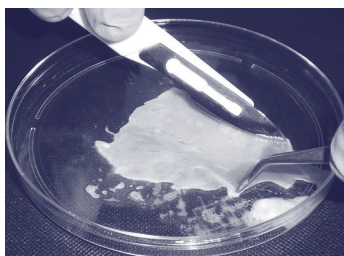

7. Subdivide skin of one mouse in two pieces.
8. Prepare enzyme mix by adding 100 µL of Enzyme G, ACF to 3.9 of mL PBS buffer in a 50 mm petri dish. Mix well.

9. Place tissue pieces on top of the enzyme mix with the dermal side facing downwards.

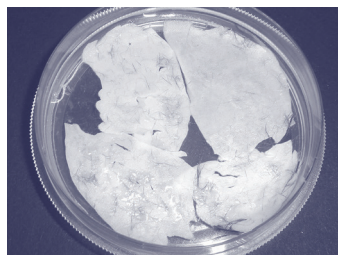


10. Incubate at 4 °C for 16 hours.


▲ **Note:** Do not extend incubation time!

2.2.2 Dissociation of epidermis (day 2)

1. Prewarm the water bath to 37 °C.
2. Peel off the epidermis from the dermis using curved tweezers.

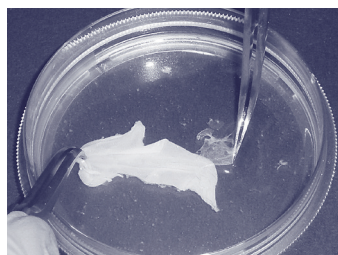

3. Transfer the epidermis into a new 50 mm petri dish containing 4 mL of 1× Buffer S and cut into smaller pieces.

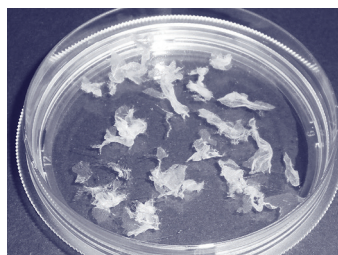

4. For automatic dissociation using the gentleMACS Dissociator proceed with chapter 2.2.3.

2.2.3 Automated dissociation using the gentleMACS Dissociator

1. Prepare enzyme mix by adding 3.9 mL of 1× Buffer S, 100 µL of Enzyme P, ACF, and 5 µL of Enzyme A, ACF into a gentleMACS C Tube.
2. Transfer epidermis pieces into the gentleMACS C Tube containing the enzyme mix. The sample material must not stick to the wall of the tube.
3. Incubate for 20 minutes at 37 °C in a water bath.

▲ **Note:** Ensure that the sample material is located in the enzyme mix during the incubation time.

4. Stop enzymatic reaction by adding 4 mL of PB buffer.

5. Tightly close C Tube and attach it upside down onto the sleeve of the gentleMACS Dissociator.

▲ **Note:** Close C Tube tightly beyond the first resistance.

▲ **Note:** Ensure that the sample material is located in the enzyme mix in the area of the rotor/stator before starting the gentleMACS Program.

6. Run the gentleMACS Program B.
7. After termination of the program, detach C Tube from the gentleMACS Dissociator.
8. Apply sample to a MACS SmartStrainer (70 µm) placed on a 50 mL tube.
9. Wash the MACS SmartStrainer (70 µm) with 10 mL of PB buffer.
10. Discard the MACS SmartStrainer (70 µm) and centrifuge sample at 300×g for 10 minutes at room temperature.
11. Aspirate supernatant completely and resuspend cells with PB buffer or an adequate buffer to the required volume for further applications.
 ▲ **Note:** Do not vortex.
 ▲ **Note:** If cell clumps occur after the washing step, add another 2.5 µL of Enzyme A, ACF per mL of cell suspension. Mix gently and incubate for 5 minutes at 37 °C in a water bath. Centrifuge at 300×g for 10 minutes and repeat step 11.
12. Process cells immediately for further applications.

Refer to www.miltenyibiotec.com for all data sheets and protocols. Miltenyi Biotec provides technical support worldwide. Visit www.miltenyibiotec.com for local Miltenyi Biotec Technical Support contact information.

Legal notices

Limited product warranty

Miltenyi Biotec B.V. & Co. KG and/or its affiliate(s) warrant this product to be free from material defects in workmanship and materials and to conform substantially with Miltenyi Biotec's published specifications for the product at the time of order, under normal use and conditions in accordance with its applicable documentation, for a period beginning on the date of delivery of the product by Miltenyi Biotec or its authorized distributor and ending on the expiration date of the product's applicable shelf life stated on the product label, packaging or documentation (as applicable) or, in the absence thereof, ONE (1) YEAR from date of delivery ("Product Warranty"). Miltenyi Biotec's Product Warranty is provided subject to the warranty terms as set forth in Miltenyi Biotec's General Terms and Conditions for the Sale of Products and Services available on Miltenyi Biotec's website at www.miltenyibiotec.com, as in effect at the time of order ("Product Warranty"). Additional terms may apply. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING IF A PRODUCT IS SUITABLE FOR CUSTOMER'S PARTICULAR PURPOSE AND APPLICATION METHODS.

Technical information

The technical information, data, protocols, and other statements provided by Miltenyi Biotec in this document are based on information, tests, or experience which Miltenyi Biotec believes to be reliable, but the accuracy or completeness of such information is not guaranteed. Such technical information and data are intended for persons with knowledge and technical skills sufficient to assess and apply their own informed judgment to the information. Miltenyi Biotec shall not be liable for any technical or editorial errors or omissions contained herein.

All information and specifications are subject to change without prior notice. Please contact Miltenyi Biotec Technical Support or visit www.miltenyibiotec.com for the most up-to-date information on Miltenyi Biotec products.

Licenses

This product and/or its use may be covered by one or more pending or issued patents and/or may have certain limitations. Certain uses may be excluded by separate terms and conditions. Please contact your local Miltenyi Biotec representative or visit Miltenyi Biotec's website at www.miltenyibiotec.com for more information.

The purchase of this product conveys to the customer the non-transferable right to use the purchased amount of the product in research conducted by the customer (whether the customer is an academic or for-profit entity). This product may not be further sold. Additional terms and conditions (including the terms of a Limited Use Label License) may apply.

CUSTOMER'S USE OF THIS PRODUCT MAY REQUIRE ADDITIONAL LICENSES DEPENDING ON THE SPECIFIC APPLICATION. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING FOR ITSELF WHETHER IT HAS ALL APPROPRIATE LICENSES IN PLACE. Miltenyi Biotec provides no warranty that customer's use of this product does not and will not infringe intellectual property rights owned by a third party. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

Trademarks

gentleMACS, MACS, and the Miltenyi Biotec logo are registered trademarks or trademarks of Miltenyi Biotec B.V. & Co. KG and/or its affiliates in various countries worldwide.

Copyright © 2023 Miltenyi Biotec and/or its affiliates. All rights reserved.