

## Contents

1. Description
  - 1.1 Principle of the Epidermis Dissociation Kit
  - 1.2 Background information
  - 1.3 Applications
  - 1.4 Reagent and instrument requirements
2. Protocols
  - 2.1 Reagent and instrument preparation
  - 2.2 Epidermis dissociation protocol
    - 2.2.1 Separation of epidermis from dermis (day 1)
    - 2.2.2 Automated dissociation of epidermis using the gentleMACS™ Dissociator (day 2)
3. Appendix

## 1. Description

<b>Components</b>	3 vials, containing <b>1 vial of Enzyme G (lyophilized powder)</b> <b>2.5 mL of Enzyme P</b> <b>1 vial of Enzyme A (lyophilized powder)</b>
<b>Size</b>	For 100 digestions. The specified number of digestions is valid for 4 mm biopsies following the protocol in chapter 2.2.
<b>Storage</b>	Upon arrival immediately store Enzyme P in aliquots at -20 °C. Store all other components at 2–8 °C upon arrival. Reconstitute all components before the date indicated on the box label. For information about reconstitution and storage after reconstitution of the lyophilized components refer to chapter 2.1.

### 1.1 Principle of the Epidermis Dissociation Kit

Human epidermal tissue can be dissociated to single-cell suspensions by combining mechanical dissociation with enzymatic degradation of the extracellular adhesion proteins, which maintain the structural integrity of tissues.

In a first step, the epidermal cell layer is removed from the dermal cell layer after enzymatic treatment over night at 4 °C. The epidermal tissue is then further digested enzymatically and dissociated into a single-cell suspension by using the gentleMACS™ Dissociators. Cells should be processed immediately for downstream applications, such as cell separation, cellular or molecular analyses.

### 1.2 Background information

The Epidermis Dissociation Kit, human enables the gentle and efficient generation of single-cell suspensions from human epidermal tissue. The kit has been particularly developed for the isolation of keratinocytes from diverse human skin biopsies. Furthermore, dissociated cells can be subsequently cultured or isolated using MACS® Technology.

### 1.3 Applications

- Dissociation of human skin tissue for the cultivation of keratinocytes.
- Phenotyping or enumeration of human epidermal cell populations by flow cytometry.

### 1.4 Reagent and instrument requirements

- RPMI 1640 (# 130-091-440)
- PBS: phosphate-buffered saline pH 7.2
- PB buffer: Prepare a solution containing phosphate-buffered saline (PBS), pH 7.2, 0.5% bovine serum albumin (BSA) by diluting MACS BSA Stock Solution (# 130-091-376) 1:20 with PBS. Keep buffer cold (2–8 °C).
- Pre-Separation Filter, 70 µm (# 130-095-823)
- MACSmix™ Tube Rotator (# 130-090-753) in combination with an incubator at 4 °C
- gentleMACS Dissociator (# 130-093-235) or gentleMACS Octo Dissociator (# 130-095-937)
- gentleMACS C Tubes (# 130-093-237, # 130-096-334)
- (Optional) MACS Tissue Storage Solution (# 130-100-008)
- (Optional) Tool for taking punch biopsies (e.g., Biopsy Punch)
- (Optional) ART® 1000 REACH™ pipet tips (Molecular BioProducts, Inc.) for removal of dissociated material from the closed C Tubes.

## 2. Protocols

▲ For details on the use of the gentleMACS Dissociators, refer to the gentleMACS Dissociator user manuals.

▲ For cell culture experiments subsequent to tissue dissociation, all steps should be performed under sterile conditions.

▲ The protocol has been optimized for the digestion of adult human skin from breast or abdominal reduction surgery.

▲ Up to five punch biopsies (4 mm each) can be used per digestion. When working with bigger punch biopsies, cut the biopsy into pieces with a maximum diameter of 4 mm.

### 2.1 Reagent and instrument preparation

▲ Prepare Enzyme G by reconstitution of the lyophilized powder in the vial with 3 mL of sterile, distilled water. Do not vortex. Produce aliquots of appropriate volume. Store aliquots at  $-20^{\circ}\text{C}$ . Avoid repeated freeze-thaw-cycles.

▲ Prepare aliquots of appropriate volume of Enzyme P. Store aliquots at  $-20^{\circ}\text{C}$ . Avoid repeated freeze-thaw-cycles.

▲ Prepare Enzyme A by reconstitution of the lyophilized powder in the vial with 1 mL RPMI 1640. Do not vortex. Produce aliquots of appropriate volume. Store aliquots at  $-20^{\circ}\text{C}$ . Avoid repeated freeze-thaw-cycles.

### 2.2 Epidermis dissociation protocol

#### 2.2.1 Separation of epidermis from dermis (day 1)

1. Wash human skin tissue in an appropriate buffer or cell culture medium, e.g., MACS Tissue Storage Solution.
2. Remove subcutaneous fat using scissors. If the diameter of the skin sample exceeds 4 mm in diameter, take one or more 4 mm diameter punches by rotating down the tool (e.g., Biopsy Punch) through epidermis and dermis. Store punches in an appropriate buffer or cell culture medium, e.g., MACS Tissue Storage Solution, until needed.
3. Transfer 1 mL of RPMI 1640 and 25  $\mu\text{L}$  of Enzyme G to into a 2 mL tube and mix carefully. Keep on ice.
4. Transfer up to five samples of skin tissue (4 mm diameter) into the 2 mL tube.
5. Incubate sample for 14–18 hours at  $4^{\circ}\text{C}$  using the MACSmix Tube Rotator (12 rpm).

#### 2.2.2 Automated dissociation of epidermis using the gentleMACS™ Dissociator (day 2)

1. Prewarm the water bath to  $37^{\circ}\text{C}$ .
2. Take the biopsy out of the 2 mL tube and peel off the epidermis from the dermis using curved tweezers.
3. (Optional) Transfer the epidermis without dermis back to the 2 mL tube and incubate for additional 2 hours at  $37^{\circ}\text{C}$ .  
▲ **Note:** This step will increase yield of Langerhans cells.
4. Prepare enzyme mix by adding 1 mL of RPMI 1640, 25  $\mu\text{L}$  of Enzyme P, and 5  $\mu\text{L}$  of Enzyme A into a new 2 mL tube. Keep on ice.  
▲ **Note:** Do not mix Enzyme P and Enzyme A directly.
5. Transfer the epidermis into the 2 mL tube containing the enzyme mix.

6. Incubate for 60 minutes at  $37^{\circ}\text{C}$  in a water bath.

▲ **Note:** It has to be ensured that the sample material is located in the enzyme mix during the incubation time.

7. Lay the epidermis directly on the rotator/stator of the gentleMACS™ C Tube. Transfer the remaining liquid of the 2 mL tube in the C Tube.

▲ **Note:** It has to be ensured that the sample material is located in the enzyme mix in the area of the rotor/stator before starting the gentleMACS Program as the epidermis might easily stick to the tube wall and won't be dissociated.

8. Stop enzymatic reaction by adding 1 mL of cold PB buffer.
9. Tightly close C Tube and attach it upside down onto the sleeve of the gentleMACS Dissociator.

▲ **Note:** Close C Tube tightly beyond the first resistance.

10. Run the gentleMACS Program B.

11. After termination of the program, detach C Tube from the gentleMACS Dissociator.

12. Perform a short centrifugation step to collect the sample material at the tube bottom.

13. Resuspend sample by pipetting up and down and apply the cell suspension to a Pre-Separation Filter, 70  $\mu\text{m}$  placed on a 15 mL tube.

▲ **Note:** Dissociated tissue can be removed from the closed C Tube by pipetting through the septum-sealed opening in the center of the cap of the C Tube. Use ART 1000 REACH 1000  $\mu\text{L}$  pipette tips.

14. Wash the Pre-Separation Filter, 70  $\mu\text{m}$ , with 1 mL of cold PB Buffer.

▲ **Note:** (Optional) To collect remaining cells in the C Tube add buffer first to the C Tube and then on top of the filter.

15. Discard the Pre-Separation Filter, 70  $\mu\text{m}$ , and centrifuge sample at  $300\times g$  for 10 minutes at room temperature. Aspirate supernatant completely

16. Resuspend cells by pipetting up and down with PB buffer or an appropriate buffer to the required volume for further applications.

▲ **Note:** If cell clumps occur after the washing step, add another 10  $\mu\text{L}$  of Enzyme A per mL of cell suspension, mix gently, incubate for 5 minutes at  $37^{\circ}\text{C}$  in a water bath, and centrifuge at  $300\times g$  for 10 minutes. Aspirate supernatant completely and repeat step 16.

17. Process cells immediately for further applications.

### 3. Appendix

#### Target cell yield per sample

Typical target cell yields per 4 mm punch biopsy of adult abdominal skin are about  $8 \times 10^4$  total cells (mostly keratinocytes) and  $3 \times 10^3$  Langerhans cells.

#### Typical staining

Epitopes which are intact after the dissociation procedure are, e.g., CD29 and CD49f (both positive for keratinocytes), CD45, CD207, CD1a, and HLA-DR (all positive for Langerhans cells).

Epitopes which are degraded after the dissociation procedure are, for example, CD117 (marker for melanocytes).

Refer to [www.miltenyibiotec.com](http://www.miltenyibiotec.com) for all data sheets and protocols. Miltenyi Biotec provides technical support worldwide. Visit [www.miltenyibiotec.com](http://www.miltenyibiotec.com) for local Miltenyi Biotec Technical Support contact information.

### Legal notices

#### Limited product warranty

Miltenyi Biotec B.V. & Co. KG and/or its affiliate(s) warrant this product to be free from material defects in workmanship and materials and to conform substantially with Miltenyi Biotec's published specifications for the product at the time of order, under normal use and conditions in accordance with its applicable documentation, for a period beginning on the date of delivery of the product by Miltenyi Biotec or its authorized distributor and ending on the expiration date of the product's applicable shelf life stated on the product label, packaging or documentation (as applicable) or, in the absence thereof, ONE (1) YEAR from date of delivery ("Product Warranty"). Miltenyi Biotec's Product Warranty is provided subject to the warranty terms as set forth in Miltenyi Biotec's General Terms and Conditions for the Sale of Products and Services available on Miltenyi Biotec's website at [www.miltenyibiotec.com](http://www.miltenyibiotec.com), as in effect at the time of order ("Product Warranty"). Additional terms may apply. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING IF A PRODUCT IS SUITABLE FOR CUSTOMER'S PARTICULAR PURPOSE AND APPLICATION METHODS.

#### Technical information

The technical information, data, protocols, and other statements provided by Miltenyi Biotec in this document are based on information, tests, or experience which Miltenyi Biotec believes to be reliable, but the accuracy or completeness of such information is not guaranteed. Such technical information and data are intended for persons with knowledge and technical skills sufficient to assess and apply their own informed judgment to the information. Miltenyi Biotec shall not be liable for any technical or editorial errors or omissions contained herein.

All information and specifications are subject to change without prior notice. Please contact Miltenyi Biotec Technical Support or visit [www.miltenyibiotec.com](http://www.miltenyibiotec.com) for the most up-to-date information on Miltenyi Biotec products.

#### Licenses

This product and/or its use may be covered by one or more pending or issued patents and/or may have certain limitations. Certain uses may be excluded by separate terms and conditions. Please contact your local Miltenyi Biotec representative or visit Miltenyi Biotec's website at [www.miltenyibiotec.com](http://www.miltenyibiotec.com) for more information.

The purchase of this product conveys to the customer the non-transferable right to use the purchased amount of the product in research conducted by the customer (whether the customer is an academic or for-profit entity). This product may not be further sold. Additional terms and conditions (including the terms of a Limited Use Label License) may apply.

CUSTOMER'S USE OF THIS PRODUCT MAY REQUIRE ADDITIONAL LICENSES DEPENDING ON THE SPECIFIC APPLICATION. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING FOR ITSELF WHETHER IT HAS ALL APPROPRIATE LICENSES IN PLACE. Miltenyi Biotec provides no warranty that customer's use of this product does not and will not infringe intellectual property rights owned by a third party. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

#### Trademarks

gentleMACS, MACS, MACSmix, and the Miltenyi Biotec logo are registered trademarks or trademarks of Miltenyi Biotec and/or its affiliates in various countries worldwide. All other trademarks mentioned in this publication are the property of their respective owners and are used for identification purposes only.

Copyright © 2021 Miltenyi Biotec and/or its affiliates. All rights reserved.