

Contents

1. Description
 - 1.1 Principle of the MACS® Separation
 - 1.2 Background information
 - 1.3 Applications
 - 1.4 Reagent and instrument requirements
2. Protocol
 - 2.1 Sample preparation
 - 2.2 Magnetic labeling of non-endothelial cells
 - 2.3 Magnetic separation: Depletion of non-endothelial cells
 - 2.4 Magnetic labeling of cardiac endothelial cells
 - 2.5 Magnetic separation: Positive selection of neonatal cardiac endothelial cells
3. Example of a separation using the Neonatal Cardiac Endothelial Cell Isolation Kit

Warnings

Reagents contain sodium azide. Under acidic conditions sodium azide yields hydrazoic acid, which is extremely toxic. Azide compounds should be diluted with running water before discarding. These precautions are recommended to avoid deposits in plumbing where explosive conditions may develop.

1. Description

This product is for research use only.

Components	1 mL Non-Cardiac Endothelial Cell Depletion Cocktail, mouse 1 mL Cardiac Endothelial Cell Isolation Cocktail, mouse
Capacity	For 2.5×10^8 total cells, up to 50 separations.
Product format	All components are supplied in buffer containing stabilizer and 0.05% sodium azide.
Storage	Store protected from light at 2–8 °C. Do not freeze. The expiration date is indicated on the vial label.

1.1 Principle of the MACS® Separation

The isolation of cardiac endothelial cells is performed in a two-step procedure. First, the non-endothelial cells are magnetically labeled with Non-Cardiac Endothelial Cell Depletion Cocktail. The labeled cells are subsequently depleted by separation over a MACS® Column, which is placed in the magnetic field of a MACS Separator. In the second step, the cardiac endothelial cells are labeled with Cardiac Endothelial Cell Isolation Cocktail and isolated by positive selection from the non-endothelial cell-depleted fraction by separation over a MACS Column, which is placed in the magnetic field of a MACS Separator.

After removing the column from the magnetic field, the magnetically retained neonatal cardiac endothelial cells can be eluted as the positively selected cell fraction.

Neonatal mouse hearts (P0–P3): Depletion of non-endothelial cells

1. Magnetic labeling of non-endothelial cells with Non-Cardiac Endothelial Cell Depletion Cocktail.
2. Magnetic separation using an LD Column or an autoMACS Column (program "Deplete05").

Pre-enriched cardiac endothelial cells (flow-through fraction): Positive selection of neonatal cardiac endothelial cells

1. Magnetic labeling of cardiac endothelial cells with Cardiac Endothelial Cell Isolation Cocktail.
2. Magnetic separation using an MS Column or an autoMACS Column (program "Possel").

Neonatal cardiac endothelial cells

1.2 Background information

Besides cardiomyocytes and fibroblasts, endothelial cells represent the third most frequent cell population in the neonatal mouse heart. Endothelial cells form the inner layer of blood vessels and play a key role in vascular development, homeostasis and remodeling. The Neonatal Cardiac Endothelial Cell Isolation Kit, mouse has been designed for enrichment of cardiac endothelial cells from dissociated neonatal mouse hearts. The isolation of cardiac endothelial cells has successfully been tested using neonatal mouse hearts from postnatal day 0 to day 3 (P0-P3). For optimal use, the Neonatal Cardiac Endothelial Cell Isolation Kit should be used in combination with the Neonatal Heart Dissociation Kit, mouse and rat (# 130-098-373).

1.3 Applications

- Enrichment of cardiac endothelial cells from neonatal (P0-P3) mouse hearts.
- Culture and expansion or direct use of enriched cardiac endothelial cells for biochemical, physiological and pharmacological studies.

1.4 Reagent and instrument requirements

- Buffer: Prepare a solution containing phosphate-buffered saline (PBS), pH 7.2, 0.5% bovine serum albumin (BSA). Keep buffer cold (2–8 °C). Degas buffer before use, as air bubbles could block the column.

▲ **Note:** Do not use ready-to-use buffers, e.g., autoMACS Running Buffer or autoMACS Rinsing Buffer.

▲ **Note:** Addition of DNase (230 U/mL) to the buffer is optional, but may improve plating of endothelial cells after cell separation.

▲ **Note:** BSA can be replaced by other proteins such as mouse serum albumin, mouse serum, or fetal bovine serum (FBS). Buffers or media containing Ca²⁺ or Mg²⁺ are not recommended for use.

- MACS Columns and MACS Separators: For optimal purity and recovery the use of an LD Column for depletion of non-cardiac endothelial cells is strongly recommended. For the subsequent positive selection of neonatal cardiac endothelial cells the use of an MS Column is strongly recommended. Positive selection and depletion can also be performed by using the autoMACS Pro or the autoMACS Separator if the use of the autoMACS Running Buffer or autoMACS Rinsing Solution is avoided. Please use buffers for separation as indicated in this section.

Column	Max. number of labeled cells	Max. number of total cells	Separator
Depletion			
LD	10 ⁸	5×10 ⁸	MidiMACS, QuadroMACS, VarioMACS, SuperMACS II
Positive selection			
MS	10 ⁷	2×10 ⁸	MiniMACS, OctoMACS, VarioMACS, SuperMACS II
Positive selection or depletion			
autoMACS	2×10 ⁸	4×10 ⁹	autoMACS Pro

▲ **Note:** Column adapters are required to insert certain columns into the VarioMACS™ or SuperMACS™ II Separators. For details refer to the respective MACS Separator data sheet.

- (Optional) Fluorochrome-conjugated antibodies for flow cytometric analysis, e.g., CD90.2-APC (# 130-091-790) or CD31-PE (# 130-097-422). For more information about antibodies refer to www.miltenyibiotec.com/antibodies.
- Neonatal Heart Dissociation Kit (# 130-098-373) for the generation of single-cell suspension from neonatal heart tissue.
- Red Blood Cell Lysis Solution (10×) (130-094-183).
- gentleMACS Dissociator (# 130-093-235), gentleMACS Octo Dissociator (# 130-095-937) or gentleMACS Octo Dissociator with Heaters (# 130-096-427), and gentleMACS C Tubes (# 130-093-237, # 130-096-334).
- (Optional) Propidium Iodide Solution (# 130-093-233) or 7-AAD for flow cytometric exclusion of dead cells.
- (Optional) Pre-Separation Filters (70 μm, # 130-095-823) to remove cell clumps.

2. Protocol

2.1 Sample preparation

For preparation of single-cell suspensions from neonatal hearts use the Neonatal Heart Dissociation Kit, mouse and rat (# 130-098-373) in combination with the gentleMACS Dissociator (# 130-092-235), the gentleMACS Octo Dissociator (# 130-095-937) or gentleMACS Octo Dissociator with Heaters (# 130-096-427).

For efficient plating and culture of isolated cardiac endothelial cells it is strongly recommended to use fibronectin-coated cell culture dishes. Coat cell culture dishes with fibronectin (100 μg/mL) at least for 2 hours in the incubator. Before use aspirate the coating solution and add the cell suspension immediately.

For optimal performance it is recommended to perform a red blood cell lysis before magnetic labeling and separation by using Red Blood Cell Lysis Solution (10×) (# 130-094-183). For details, e.g., regarding optimal incubation time, refer to the Neonatal Heart Dissociation Kit, mouse and rat data sheet.



2.2 Magnetic labeling of non-endothelial cells

▲ Work fast, keep cells cold, and use pre-cooled solutions. This will prevent capping of antibodies on the cell surface and non-specific cell labeling.

▲ Volumes for magnetic labeling given below are for up to 5×10⁶ total cells. When working with fewer than 5×10⁶ cells, use the same volumes as indicated. When working with higher cell numbers, scale up all reagent volumes and total volumes accordingly (e.g. for 10⁷ total cells, use twice the volume of all indicated reagent volumes and total volumes).

▲ For optimal performance it is important to obtain a single-cell suspension before magnetic labeling. Pass cells through 70 μm nylon mesh (Pre-Separation Filters, 70 μm, # 130-095-823) to remove cell clumps which may clog the column. Moisten filter with buffer before use.

▲ The recommended incubation temperature is 2–8 °C. Higher temperatures and/or longer incubation times may lead to non-specific cell labeling. Working on ice may require increased incubation times.

1. Determine cell number.
2. Centrifuge cell suspension at 300×g for 5 minutes. Aspirate supernatant completely.
3. Resuspend cell pellet in 80 μL of buffer per 5×10⁶ total cells.
4. Add 20 μL of Non-Cardiac Endothelial Cell Depletion Cocktail per 5×10⁶ total cells.
5. Mix well and incubate for 15 minutes in the refrigerator (2–8 °C).
6. Adjust volume to 500 μL using buffer for 5×10⁶ total cells. Do not centrifuge.
7. Proceed to magnetic separation (2.3).



2.3 Magnetic separation: Depletion of non-endothelial cells

▲ Choose an LD Column and Separator. For details refer to the table in section 1.4.

▲ Always wait until the column reservoir is empty before proceeding to the next step.

1. Place LD Column in the magnetic field of a suitable MACS Separator. For details see LD Column data sheet.
2. Prepare column by rinsing with 2 mL of buffer.
3. Apply cell suspension onto the column.
4. Collect unlabeled cells that pass through and wash column with 2×1 mL of buffer. Collect total effluent; this is the unlabeled pre-enriched endothelial cell fraction.
5. (Optional) Remove column from the separator and place it on a suitable collection tube. Pipette 3 mL of buffer onto the column. Immediately flush out the magnetically labeled non-endothelial cells by firmly pushing the plunger into the column
6. Proceed to 2.4 for the labeling of neonatal cardiac endothelial cells.

Depletion with the autoMACS® Pro Separator

▲ Refer to the respective user manual for instructions on how to use the autoMACS® Pro Separator.

▲ Buffers used for operating the autoMACS Pro Separator should have a temperature of ≥ 10 °C.

1. Prepare and prime the instrument.
2. Apply tube containing the sample and provide tubes for collecting the labeled and unlabeled cell fractions. Place sample tube in row A of the tube rack and the fraction collection tubes in rows B and C.
3. For a standard separation choose the following program:
Depletion: Deplete05
Collect negative fraction in row B of the tube rack. This fraction represents the pre-enriched cardiac endothelial cells.
4. Proceed to 2.4 for the labeling of neonatal cardiac endothelial cells.



2.4 Magnetic labeling of cardiac endothelial cells

▲ Volumes for magnetic labeling given below are for an initial starting cell number of up to 5×10^6 total cells. For higher initial cell numbers, scale up all volumes accordingly.

1. Centrifuge cell suspension at $300 \times g$ for 10 minutes. Aspirate supernatant completely.
2. Resuspend cell pellet in 80 μL of buffer.
3. Add 20 μL of Cardiac Endothelial Cell Isolation Cocktail.
4. Mix well and incubate for 15 minutes in the refrigerator ($2-8$ °C).
5. Wash cells by adding 1–2 mL of buffer per 5×10^6 initial cells and centrifuge at $300 \times g$ for 5 minutes. Aspirate supernatant completely.
6. Resuspend up to 5×10^6 cells in 500 μL of buffer.
▲ **Note:** For higher cell numbers, scale up buffer volume accordingly.
7. Proceed to magnetic separation (2.5).



2.5 Magnetic separation: Positive selection of neonatal cardiac endothelial cells

Positive selection with MS Columns



1. Place MS Column in the magnetic field of a suitable MACS Separator. For details refer to MS Column data sheet.
2. Prepare column by rinsing with 500 μL of buffer.
3. Apply cell suspension onto the column. Collect flow-through containing unlabeled cells.
4. Wash column with 3×500 μL of buffer. Collect unlabeled cells that pass through and combine with the effluent from step 3.
▲ **Note:** Perform washing steps by adding buffer aliquots as soon as the column reservoir is empty.
5. Remove column from the separator and place it on a suitable collection tube.
6. Pipette 1 mL of buffer onto the column. Immediately flush out the magnetically labeled cells by firmly pushing the plunger into the column.

Positive selection with the autoMACS® Pro Separator

1. Prepare and prime the instrument.
2. Apply tube containing the sample and provide tubes for collecting the labeled and unlabeled cell fractions. Place sample tube in row A of the tube rack and the fraction collection tubes in rows B and C.
3. For a standard separation choose the following program:
Positive selection: Possel
Collect positive fraction in row C of the tube rack. This is the enriched neonatal cardiac endothelial cell fraction.

3. Example of a separation using the Neonatal Cardiac Endothelial Cell Isolation Kit

Cardiac endothelial cells were isolated from P1 CD-1 mouse hearts by using the Neonatal Cardiac Endothelial Cell Isolation Kit, one LD Column with MidiMACS™ Separator, and one MS Column with MiniMACS™ Separator. The cells were fluorescently stained with CD90.2-APC (# 130-091-790) and CD31-PE (# 130-097-422) and analyzed by flow cytometry using the MACSQuant® Analyzer. Cell debris and dead cells were excluded from the analysis based on scatter signals and propidium iodide fluorescence.

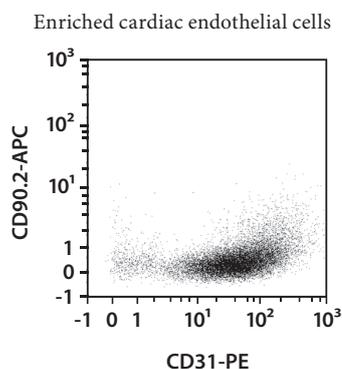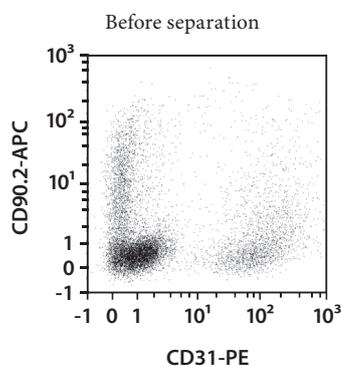

Refer to www.miltenyibiotec.com for all data sheets and protocols. Miltenyi Biotec provides technical support worldwide. Visit www.miltenyibiotec.com/local to find your nearest Miltenyi Biotec contact.

Legal notices

Limited product warranty

Miltenyi Biotec B.V. & Co. KG and/or its affiliate(s) warrant this product to be free from material defects in workmanship and materials and to conform substantially with Miltenyi Biotec's published specifications for the product at the time of order, under normal use and conditions in accordance with its applicable documentation, for a period beginning on the date of delivery of the product by Miltenyi Biotec or its authorized distributor and ending on the expiration date of the product's applicable shelf life stated on the product label, packaging or documentation (as applicable) or, in the absence thereof, ONE (1) YEAR from date of delivery ("Product Warranty"). Miltenyi Biotec's Product Warranty is provided subject to the warranty terms as set forth in Miltenyi Biotec's General Terms and Conditions for the Sale of Products and Services available on Miltenyi Biotec's website at www.miltenyibiotec.com, as in effect at the time of order ("Product Warranty"). Additional terms may apply. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING IF A PRODUCT IS SUITABLE FOR CUSTOMER'S PARTICULAR PURPOSE AND APPLICATION METHODS.

Technical information

The technical information, data, protocols, and other statements provided by Miltenyi Biotec in this document are based on information, tests, or experience which Miltenyi Biotec believes to be reliable, but the accuracy or completeness of such information is not guaranteed. Such technical information and data are intended for persons with knowledge and technical skills sufficient to assess and apply their own informed judgment to the information. Miltenyi Biotec shall not be liable for any technical or editorial errors or omissions contained herein.

All information and specifications are subject to change without prior notice. Please contact Miltenyi Biotec Technical Support or visit www.miltenyibiotec.com for the most up-to-date information on Miltenyi Biotec products.

Licenses

This product and/or its use may be covered by one or more pending or issued patents and/or may have certain limitations. Certain uses may be excluded by separate terms and conditions. Please contact your local Miltenyi Biotec representative or visit Miltenyi Biotec's website at www.miltenyibiotec.com for more information.

The purchase of this product conveys to the customer the non-transferable right to use the purchased amount of the product in research conducted by the customer (whether the customer is an academic or for-profit entity). This product may not be further sold. Additional terms and conditions (including the terms of a Limited Use Label License) may apply.

CUSTOMER'S USE OF THIS PRODUCT MAY REQUIRE ADDITIONAL LICENSES DEPENDING ON THE SPECIFIC APPLICATION. THE CUSTOMER IS SOLELY RESPONSIBLE FOR DETERMINING FOR ITSELF WHETHER IT HAS ALL APPROPRIATE LICENSES IN PLACE. Miltenyi Biotec provides no warranty that customer's use of this product does not and will not infringe intellectual property rights owned by a third party. BY USE OF THIS PRODUCT, THE CUSTOMER AGREES TO BE BOUND BY THESE TERMS.

Trademarks

autoMACS, gentleMACS, MACS, MACSQuant, MidiMACS, the Miltenyi Biotec logo, MiniMACS, OctoMACS, QuadroMACS, SuperMACS, and VarioMACS are registered trademarks or trademarks of Miltenyi Biotec and/or its affiliates in various countries worldwide.

Copyright © 2020 Miltenyi Biotec and/or its affiliates. All rights reserved.