Miltényi Biotec

## CliniMACS Prodigy® Supplementary Bag

CliniMACS Prodigy Supplementary Bag

REF 200-076-612

**Miltényi Biotec B.V. & Co. KG**

Friedrich-Ebert-Straße 68

51429 Bergisch Gladbach | Germany

☎ +49 2204 8306-3803

CE 0123

**For U.S. distributed by:**

**Miltényi Biotec Inc.**

2303 Lindbergh Street

Auburn, CA 95602 | USA

☎ 1 800 367 6227

**For U.S. CAUTION:**

Clinical Research System

Limited by Federal (or United States)

Law to Investigational Use or

under an FDA Approval

## Table of contents

|    |                                       |    |
|----|---------------------------------------|----|
| en | Instructions for use . . . . .        | 2  |
| cs | Návod k použití . . . . .             | 4  |
| da | Brugsanvisning. . . . .               | 6  |
| de | Gebrauchsanweisung . . . . .          | 8  |
| el | Οδηγίες χρήσης . . . . .              | 10 |
| es | Instrucciones de uso . . . . .        | 12 |
| fi | Käyttöohje . . . . .                  | 14 |
| fr | Notice d'utilisation . . . . .        | 16 |
| hr | Upute za upotrebu . . . . .           | 18 |
| hu | Használati útmutató . . . . .         | 20 |
| it | Istruzioni per l'uso. . . . .         | 22 |
| nl | Gebruiksaanwijzing. . . . .           | 24 |
| no | Bruksanvisning. . . . .               | 26 |
| pl | Instrukcja używania. . . . .          | 28 |
| pt | Instruções de utilização . . . . .    | 30 |
| ro | Instrucțiuni de utilizare. . . . .    | 32 |
| ru | Инструкции по использованию . . . . . | 34 |
| sk | Návod na použitie. . . . .            | 36 |
| sl | Navodila za uporabo . . . . .         | 38 |
| sv | Bruksanvisning. . . . .               | 40 |
| tr | Kullanma talimatları. . . . .         | 42 |

# Instructions for use

## Contents

**QTY** Two CliniMACS Prodigy® Supplementary Bags

## Intended purpose

The CliniMACS Prodigy Supplementary Bag is intended to reduce the risk of contact to infectious sample material in case of a leakage of a CliniMACS Prodigy Tubing Set. The CliniMACS Prodigy Supplementary Bag must only be used in combination with the CliniMACS Prodigy.

## Precautions

The instructions for use including those described in the corresponding CliniMACS Prodigy User Manuals must be followed.

All materials that have been in contact with blood or blood components must be treated and disposed as biohazardous materials. Therefore, standard hospital or institutional requirements and national legislations must be observed.

## Warnings

### **WARNING**

**Biological hazards. Depending on the biological material used, contamination or infection may lead to severe personal injury or death. The CliniMACS Prodigy Supplementary Bag might be filled with infectious sample material. Wear protective gloves, protective clothing, and safety goggles to prevent contact with skin and eyes.**

Do not use after the use-by date printed on the bag label.

Do not use if package is damaged. Risk of contamination. If the packaging is damaged, there is the risk of contamination. Before opening, inspect the packaging for damage, puncture or tears. Use the Supplementary Bag only if package is undamaged and sealed.

Do not store blood, blood fractions or cell fractions in the Supplementary Bag.

Do not re-use. The re-use of the Supplementary Bag may lead to the endangerment of the user or third party.

## Storage

$+15^{\circ}\text{C}$   $+30^{\circ}\text{C}$  The Supplementary Bag must be stored at room temperature ( $+15^{\circ}\text{C}$  to  $+30^{\circ}\text{C}$  [ $+59^{\circ}\text{F}$  to  $+86^{\circ}\text{F}$ ]).

## Further information

The CliniMACS Prodigy Supplementary Bag has not been subjected to a sterilization process.

Any serious incident that has occurred in relation to this product should be reported to Miltenyi Biotec B.V. & Co. KG – using the contact information provided – and the competent authority of the member state in which the user of this product is established.

## Handling instructions

The CliniMACS Prodigy Supplementary Bag is a pre-installed bag placed within the compartment underneath the CentriCult Unit™ (see figure 1). Fluids accidentally leaking out of the chamber into the CentriCult Unit, the Supplementary Bag must be removed from its compartment to allow for the full capacity of the bag.

It is recommended to replace the Supplementary Bag no longer than one year after it has been installed to the CliniMACS Prodigy not exceeding the use-by date. Therefore, the installation date has to be written onto the Supplementary Bag label besides the calendar symbol.

**Note:** The Supplementary Bag compartment is quite narrow. If fluid is leaking out of the chamber into the CentriCult Unit, the Supplementary Bag must be removed from its compartment to allow for the full capacity of the bag.

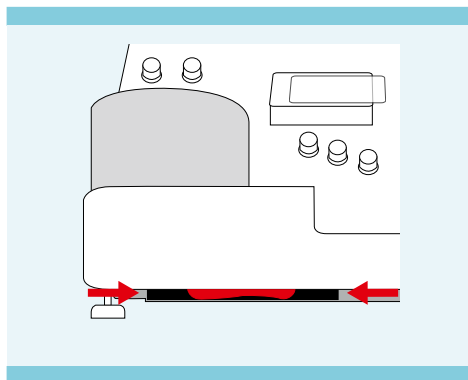
Figure 1: Supplementary Bag compartment

Let the Supplementary Bag hang down in front of the CliniMACS Prodigy until all leaking fluid is collected in the bag (see figure 2).

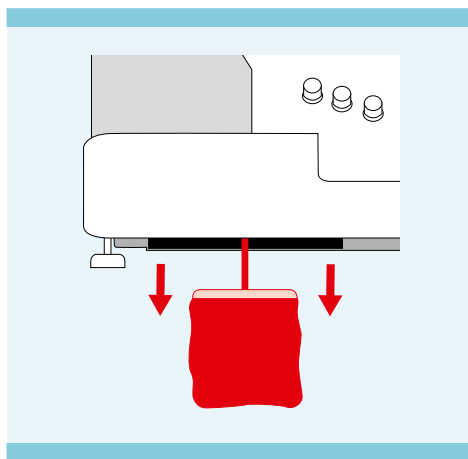
Figure 2: Collection of leaking fluid

After a leakage event the CliniMACS Prodigy must be cleaned and disinfected according to the instructions given in the CliniMACS Prodigy User Manual (Instrument).

Before starting a new process on the CliniMACS Prodigy, the filled Supplementary Bag has to be replaced by a new Supplementary Bag. One additional Supplementary Bag is delivered with the CliniMACS Prodigy. Make sure that a replacement for the Supplementary Bag is always available. It may be ordered via the Miltenyi Biotec Customer Service (REF 200-076-612).

For technical support, contact Miltenyi Biotec Technical Support.

- Clamp the tubing of the used Supplementary Bag and disconnect the bag by pulling down the connector from the tubing port under the CliniMACS Prodigy.
-  Before installation of a new Supplementary Bag, the installation date must be written on the bag label.
- Connect a new Supplementary Bag by pushing the connector of the tubing over the tubing port under the CliniMACS Prodigy.
- Make sure the tubing port is fully covered by the tubing connector (see figure 3).

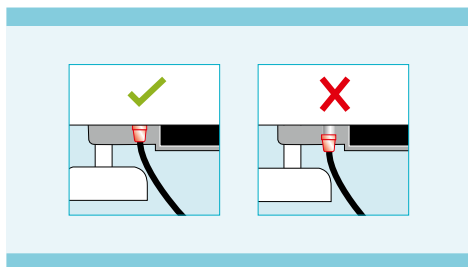
Figure 3: Correct connection of tubing connector

- Place the Supplementary Bag in the compartment underneath the CliniMACS Prodigy.
- Be careful not to obstruct the fluid pathway by kinks or bends in the tubing.

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Miltenyi Biotec Swiss AG
Miltenyi Biotec Ltd.